

NanoFrazor – A Nanolithography Tool for 2D & 3D Devices

Nils Goedecke

Jana Chabaan, ZhenMing Wu

Heidelberg Instruments Nano, Bändliweg 30, 8048 Zürich, Switzerland

Nils.goedecke@himt.ch, www.heidelberg-instruments.com

NanoFrazor systems are nanofabrication instruments based on thermal scanning probe technology. They can be used for a wide range of surface manipulation methods like nanolithography as recently summarize by Howell [1]. Core of the technology is a heatable probe tip that is used for both patterning and simultaneous inspection of complex nanostructures. The heated tip can pattern very high-resolution (< 10 nm half-pitch) nanostructures by locally evaporating resist materials. The structures are inspected by the cold tip in parallel with the patterning process, enabling stitching and markerless overlay with sub-5 nm accuracy [2]. The technique is compatible with all the common pattern transfer processes [3,4,5].

Shaping 2D materials into narrow ribbons, Hall bars etc. is often required in order to study their properties. Another challenge is formation of high-quality electrical contacts on them. Predominant fabrication process - i.e. EBL followed by etching or lift-off of metal - has its resolution limited by proximity effects, may require complex overlay procedures and typically yields poor quality non-ohmic metal contacts with high Schottky barriers and large contact resistances [6]. Here, we show that NanoFrazor lithography can be used for shaping 2D materials with very high precision (Figure 1a-b) [7] and for forming high-quality metal contact electrodes on them (Figures 1 c-d) [5]. The fabricated devices exhibit vanishing Schottky barrier heights (around 0 meV, Figure 1d).

References

- [1] Howell, S.T. *et al. Microsyst Nanoeng* **6**, 21 (2020).
- [2] Rawlings *et al.*, *ACS Nano* **9**, 6188 (2015).
- [3] Wolf *et al.*, *J. Vac. Sci. Technol. B* **33**, 02B102 (2015).
- [4] Kulmala *et al.*, *Proc. SPIE* **10584**12 (2018).
- [5] Zheng *et al.*, *Nature Electronics* **2** 17-25 (2019).
- [6] Allain *et al.*, *Nature Materials* **14**, 1195–1205 (2015).
- [7] Ryu and Knoll, *Electrical Atomic Force Microscopy for Nanoelectronics*, 143-172 (Ed. U. Celano), Springer Nature Switzerland AG (2019).

Figures

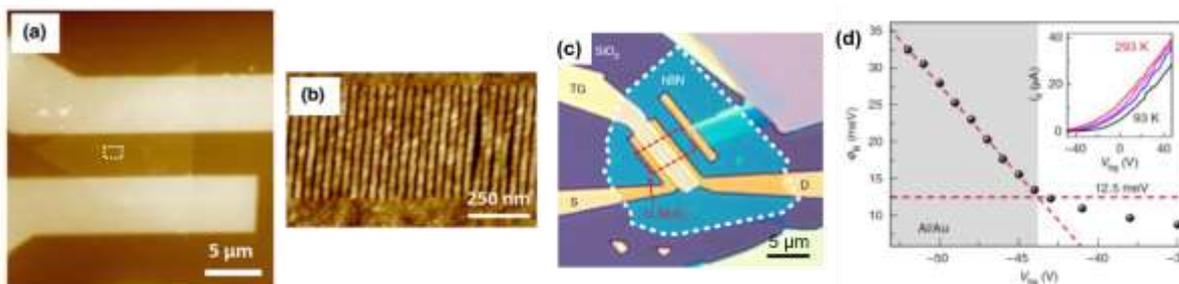


Figure 1: a) AFM image of 18-nm half-pitch 1L MoS₂ nanoribbon array patterned along the zigzag direction. (b) A close-up of the region marked with a white dashed box in (a). Figures (a) and (b) from Ref [7]. (c) Optical image of a 1L MoS₂ FET with a h-BN gate dielectric where the source, drain and top-gate electrodes have been patterned with a NanoFrazor. (d) Gate voltage dependence of Schottky barrier height of a 1L MoS₂ FET with Al/Au contacts ($V_{ds} = 2$ V). The deviation from the linear response at low V_{bg} (dashed red line) defines the flat band voltage and the SBH. Inset, corresponding temperature-dependent transfer curves ($V_{ds} = 2$ V). Figures (c) and (d) from Ref. [5].